



**ENVIROLOGIX**

For technical support call  
EnviroLogix toll free at:  
1-866-408-4597

**QuickTox™**  
*Stachybotrys/  
Aspergillus niger*  
**QuickGuide™**

*Refer to the package insert for complete instructions and interpretation*

- 1** Carefully add 15 drops of  
Extraction Buffer



- 2** Collect sample...

**For mold colonies and  
dust / particulate samples**

or

**For bulk samples such as  
drywall, ceiling tile, etc.**

*Gently swab the area until  
material is visible on swab  
head or collection device;  
insert into tube and twirl*

*Add a thumbtack-sized  
amount of material  
to the tube;  
use a pestle to grind*





- 3** Place strip into the  
extraction tube and...



Read the results!

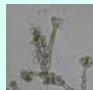

If the QuickTox Strip shows:	The result is:
no Test Line, only the Control Line	negative for <i>Stachybotrys chartarum</i> and <i>Aspergillus niger</i>
a <u>pink</u> Test Line	 positive for <i>Stachybotrys chartarum</i>
a <u>greyish-brown to greyish-black</u> Test Line	 positive for <i>Aspergillus niger</i>
a <u>brownish-pink</u> Test Line	positive for both <i>Stachybotrys chartarum</i> and <i>Aspergillus niger</i>