

QuickStix™ Kits for Soybeans
 QuickComb™ Kit for Soybean Bulk Grain
 AQ 010 BGB, AQ 014 BGB, AQ 076 TC

QuickGuide

(Refer to product inserts for detailed instructions)



Determine sample weight and grinding method. Weigh sample into an appropriate container.




Grind soybeans until finely ground. Product insert gives grind times for various methods.




Measure water using the Soybean Common Extraction™ (1:5). Add to container, cover and shake to wet entire sample.



With pipette, remove 12 mL sample from liquid on top (about 3 pipettefuls, avoid particulates), dispense into reaction cup.

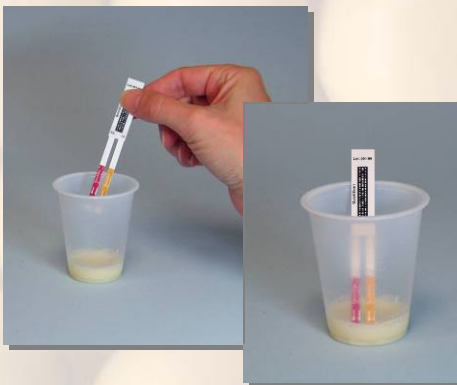

Insert comb or strip into the liquid in cup. Wait 5 minutes, then read the results—either visually or with QuickScan.

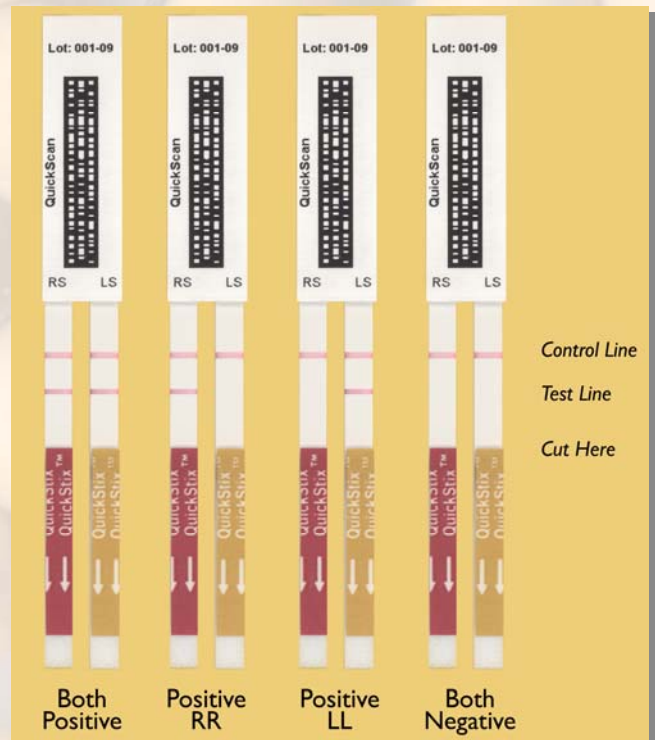

500 Riverside Industrial Parkway • Portland, ME 04103
 www.envirolgix.com • 1-866-408-4597 • 1-207-797-0300