



Sample Preparation

Prepare Solutions

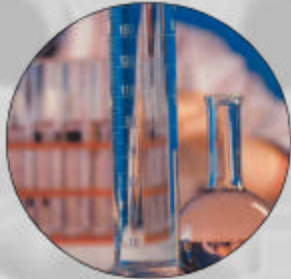

Collect & Extract Sample

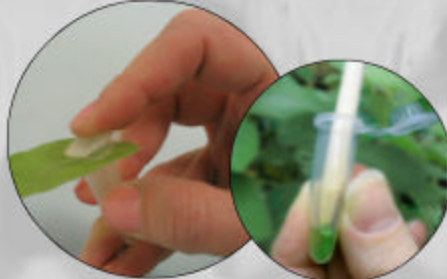


EnviroLogix


QualiPlate Kit for Soybean Rust QuickGuide

Testing Procedure


1 Add Samples, Buffer Blank,
Positive & Negative Controls

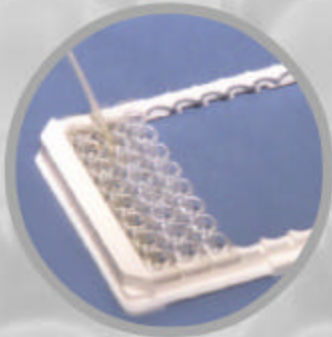


2 Mix & Incubate
for One Hour

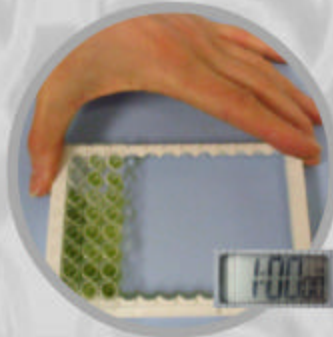


3 Wash Plate

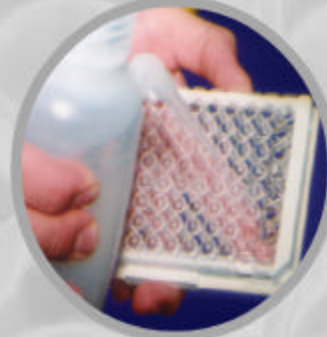


4 Add Conjugate, Mix &
Incubate for One Hour

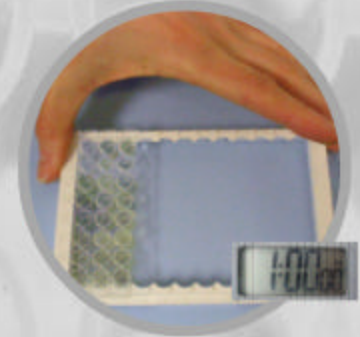


5 Wash Plate

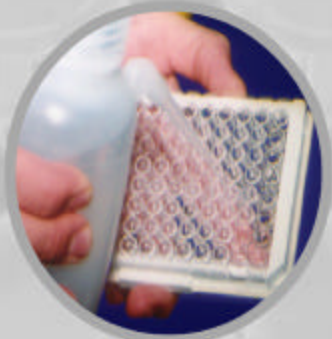

6 Add Substrate

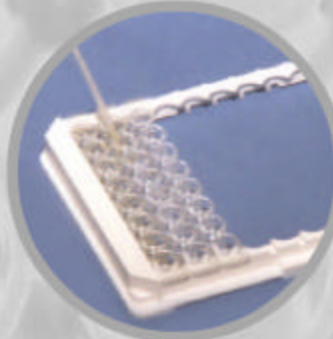

7 Mix & Incubate 20 Minutes
Read Results Visually

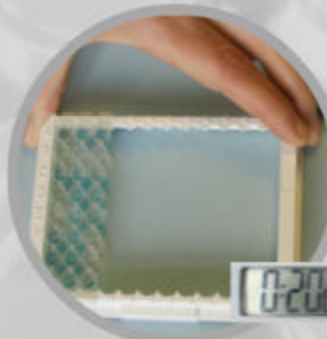

8 *OPTIONAL*~Add Stop Solution
Read Results on Reader



EnviroLogix Technical Support: Toll Free 1-866-408-4597

© EnviroLogix Inc. 2005