


QuickGuide for QuickScan

1. **Double-click the QuickScan icon** on the computer desktop -- Main Menu appears
2. After cutting tailpads off strip(s), **pull the carrier out** of the scanner to reveal the hinged compression plate (remember, strips or combs must be read immediately after reaction time)
3. Lift the plate, **place the comb or strip face-down** into the carrier, and slide it back into scanner
4. **Click “Read Test”** button on the Main Menu screen (scanner will read comb)
5. When the Results Screen appears, **enter Sample ID and/or Supplier data** where indicated (i.e. scale-ticket number or “ABC Trk 27”)
6. **Click “Save Report”** button to save the results (PDF - stored in a reports folder and accessed by a shortcut on the PC desktop)
7. **Click “Print Report”** button to create an immediate hard copy of test results, if required (“Print Receipt” prints abbreviated Report)
8. **Click “Close”** button to exit results screen and perform the next test (be sure to enter Sample ID info, or the data will not be saved)

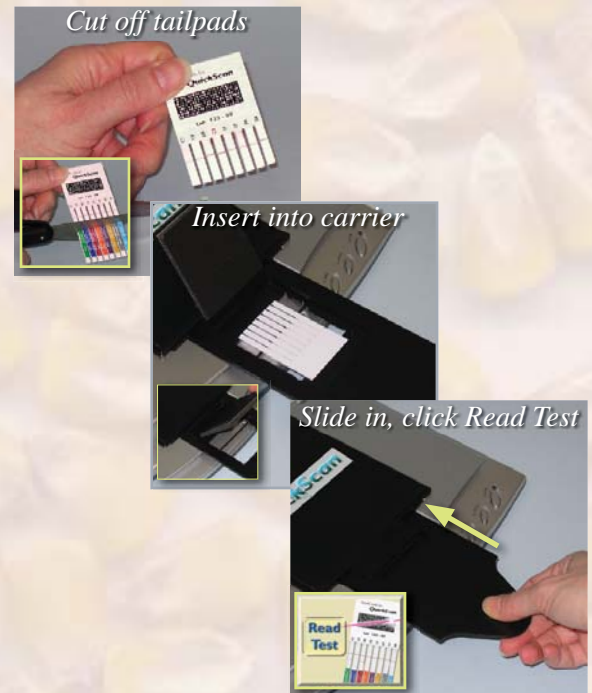

MAIN MENU



Enter Sample ID...Supplier...Comments...Action

Sample ID	Supplier	Comment 1	Comment 2	Action	Test kit	Analyte	Result(%)	TL	CL	Lot
1	Truck ABC	Farmer Bob	Corn	Very High	Accept	AQ-036 TC 13-6 C1: Cry1Ab	< 1.00			3
2						RR: Cry4E/SPS	0.94			
3						C3: Cry3Bb	1.5			
4						1F: Cry1F	2.2			
5						LP: PAT(pat)	1.8			
6						3A: Cry3A	3.7			
7						3A: mCry3A	1.5			
8						CNO Sum =	11.6			

RESULTS SCREEN