

QuickStix™ Kit for LibertyLink® Rice QuickGuide

AS-013-RB



Weigh 75 rice grains and add to blender jar; assemble with blade and place on blender base



Grind for 15-20 seconds on High; make sure all seeds are finely ground



Transfer ground rice to a weigh boat or weigh paper



Carefully pour ground rice into extraction vial



Add 1.5 times sample weight of buffer (in mL) and shake for 15 seconds



Tap solids down, place test strip in liquid portion (do not push strip into settled solids)

Wait **10 minutes**, remove strip and read result. Any clearly discernable pink Test Line is a positive result.

NOTE: Test Line intensity is dependent on GM concentration and may appear lighter than the photograph.

