

QuickStix™ Kit for Soybean Rust

QuickGuide

Sample Preparation



Tear a leaf sample



Use the QuickStix tube as a template



Leaf size is about 1 inch (2.5 cm) in diameter

Test Procedure



1. Place leaf in mesh bag and rub with coin



2. Add one pipette (5mL) of buffer



3. Massage the buffer and leaf well



4. Draw up extract with pipette




5. Fill vial with extract



6. Add a test strip to vial, wait 10 minutes

Test Interpretation

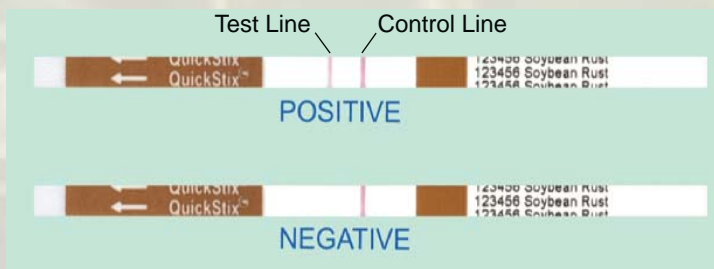

Read the result.
ANY pink Test Line is positive for SBR.

A Control Line must always be present