
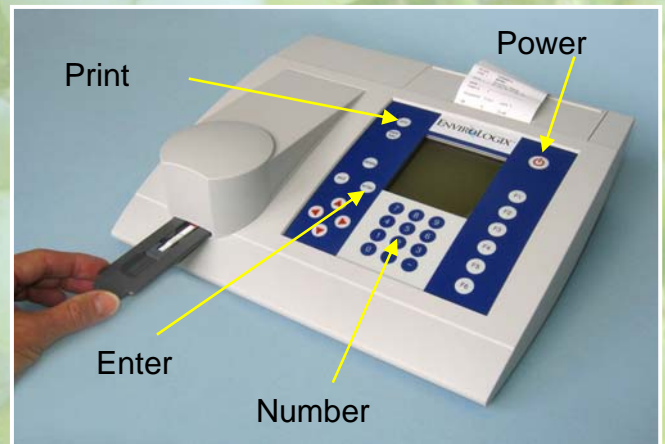


QuickStix™ Reader QuickGuide

1. **Turn on reader**
(wait 20 sec for self test/calibration)
2. Press **Enter** - (run test)
3. **Remove pad** from bottom of strip and place QuickStix into strip holder
4. **Insert strip holder** into slot and gently click into place
5. **Enter Sample ID**
(transmitted from computer if enabled);
Operator ID and Location ID are carried
OVER (change only if necessary)
6. Press **Enter**. Reader will scan barcode
and results. (if strip has no barcode, press **Enter**,
select **measurement scale** from list, press **Enter**, type in
lot #, press **Enter**)
7. Contents of the results screen will print
automatically (unless disabled, then press **Print**. See
main menu, pg.4 of manual to change setting)
and will be transmitted to your computer
(if enabled)

Test results are displayed according to Botrytis levels set by the Dewey I-W Standard (right).

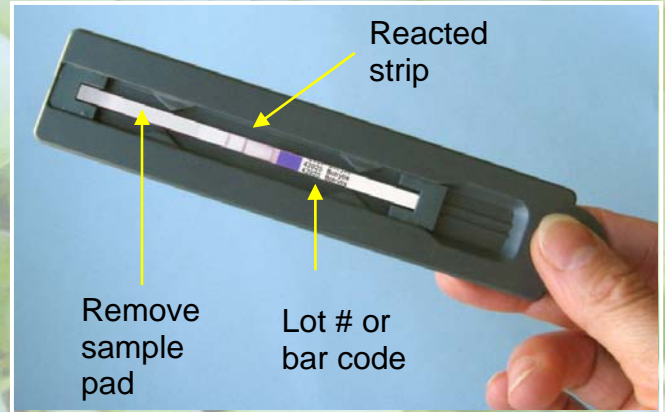

The Dewey I-W Standard

Signal Intensity	% Infection, by Incidence	% Infection, by Weight
<10	< 1.25	< 0.42
10 - 23.5	≥ 1.25 to < 2.50	≥ 0.42 to < 0.83
23.5 - 31.0	≥ 2.50 to < 5.00	≥ 0.83 to < 1.67
31.0 - 42.5	≥ 5.00 to < 10.00	≥ 1.67 to < 3.33
>42.5	≥ 10.00	≥ 3.33